



Address:
JAMBU PERSHAD & SONS
6275/22 Nicholson Road,
Ambala Cantt, Haryana,
INDIA
Pin: 133001

Email:
sales@japson.com
japsonambala@yahoo.com

Website:
www.japson.com
Phone:
+91-171-4006897

Sterility Test Apparatus

Product Image



Description

This Apparatus is made in total 316 Stainless Steel.

This apparatus is used for filtering liquid samples. The sterility tests in this system is done by the Bacteria-Retentive membranes filter method.

A 47 mm membrane filter is installed in a suitable holder. The liquid sample is drawn through this sterile membrane filter to concentrate any contaminating micro-organisms on the filter surface. After filtration is completed, sterile flushing fluid is passed into the funnels, rinsing all sample residues through the filters so that it cannot possibly inhibit growth of the retained micro-organisms. After this, the filter is sectioned and transferred to media for culturing. If growth is not visible after the incubations period, the sample passes the test.

-

Sterility Test Apparatus

Catalog No.100672

Disclaimer

The Products details given on this page are indicative in nature and JAPSON reserves the right to change them without prior notice. Buyer is also requested to re-check the specifications and other features of product at the time of order as product development is a continuous process and minor modifications may be made to design based on latest availability, process and design.